

## Automated Glucose Control in Ambr<sup>®</sup> Bioreactors Using Raman Spectroscopy

Michael Sibley<sup>1</sup>, Angus Woodhams<sup>1</sup>, Marek Hoehse<sup>2</sup>, Barney Zoro<sup>1\*</sup>

<sup>1</sup>Sartorius Royston, UK. <sup>2</sup>Sartorius Göttingen, Germany. \*barney.zoro@sartorius.com

### Aims

- Demonstrate physical integration of and automated data integration for Raman spectroscopy with Ambr<sup>®</sup> bioreactors
- Generate data to build a Raman model for glucose concentration in a typical cell culture
- Build a Raman model for glucose concentration using SIMCA 16 software
- Demonstrate real-time feedback control of glucose in culture based on the Smbr<sup>®</sup> derived SIMCA 16 Raman model

### 1. BioPAT<sup>®</sup> Spectro in Ambr<sup>®</sup> Overview

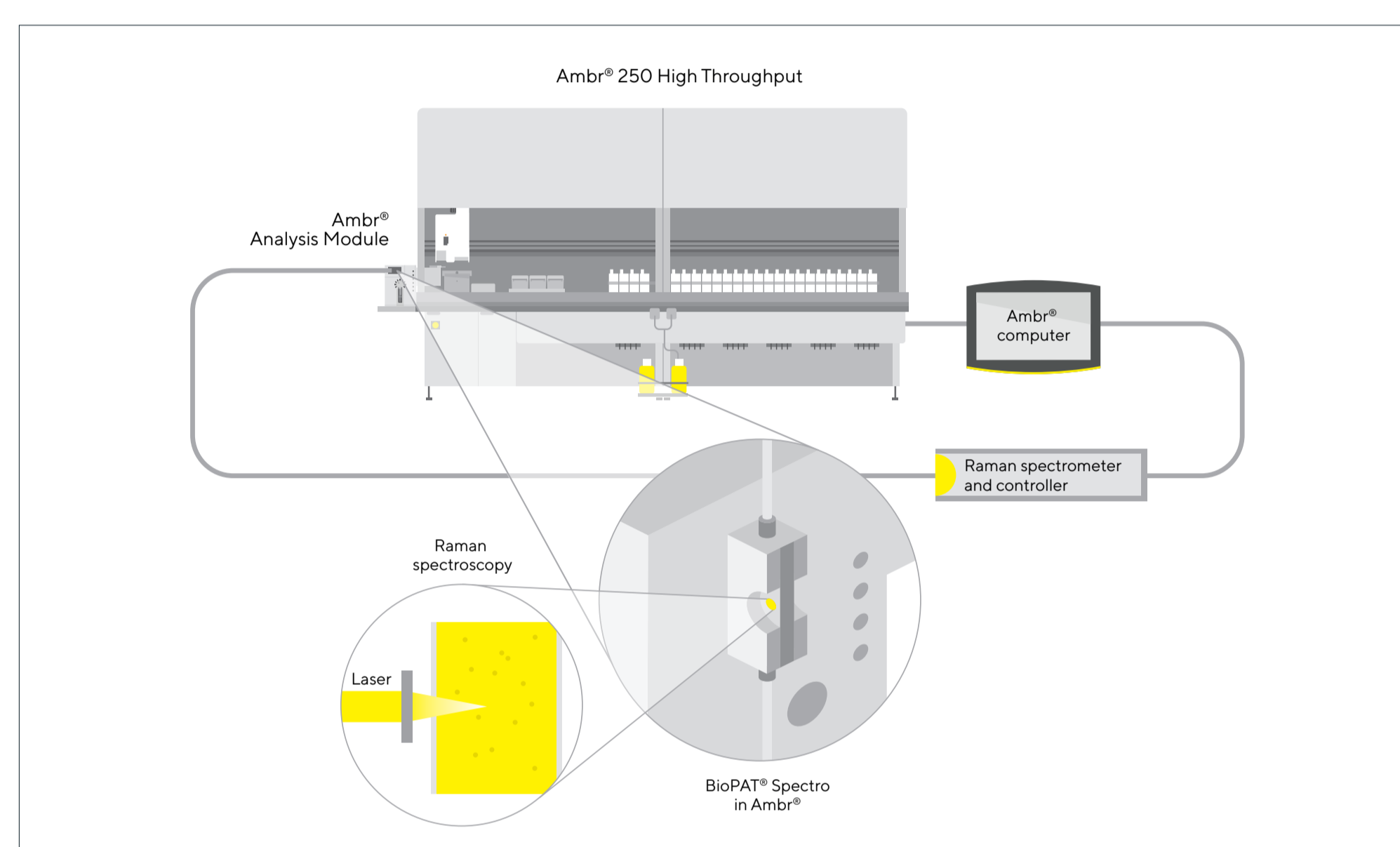


Figure 1. Schematic diagram of Ambr<sup>®</sup> Analysis Module with BioPAT<sup>®</sup> Spectro, fitted to an Ambr<sup>®</sup> 250 High Throughput bioreactor system

### 2. BioPAT<sup>®</sup> Spectro in Ambr<sup>®</sup> Key Features

- A user-replaceable Raman flow cell is integrated in the Ambr<sup>®</sup> Analysis Module, compatible with both Ambr<sup>®</sup> 15 Cell Culture and Ambr<sup>®</sup> 250 High Throughput
- Compatible with Raman Spectrometers and BioPAT<sup>®</sup> Spectro Raman probes from Kaiser Optical Systems and Tornado Spectral Systems
- Small sample volumes: 160 µL for Ambr<sup>®</sup> 15, 200 µL for Ambr<sup>®</sup> 250 High Throughput
- Fast cycle time: ~10-15 min per sample
- Samples can be automatically spiked with a concentrated analyte stock solution prior to Raman analysis
- All key data from the Ambr<sup>®</sup> process, Raman and other integrated analyzers are automatically collated by the Ambr<sup>®</sup> Software
- Identical Raman probes and flow cell optical paths support direct model transfer to Biostat STR<sup>®</sup>
- Requires compatible Ambr<sup>®</sup> Analysis Module, Win10 Ambr<sup>®</sup> Control PC, separate SIMCA 16 license for creating models
- Suitable for typical mammalian fed-batch cell densities; microbial applications are not supported


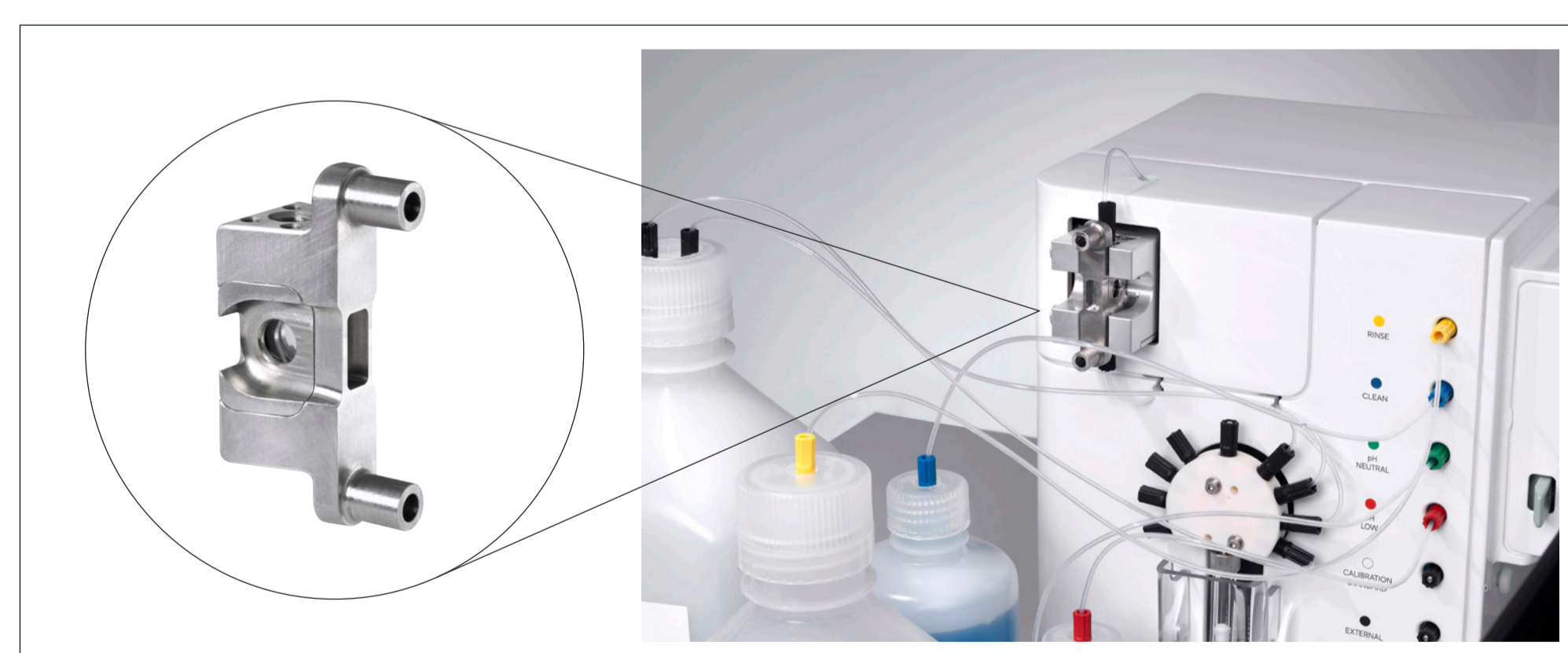


Figure 2. Ambr<sup>®</sup> Analysis Module with BioPAT<sup>®</sup> Spectro fitted to an Ambr<sup>®</sup> 15 system. Inset box: User replaceable BioPAT<sup>®</sup> Spectro flow cell.

Workflow diagram for BioPAT<sup>®</sup> Spectro in Ambr<sup>®</sup>

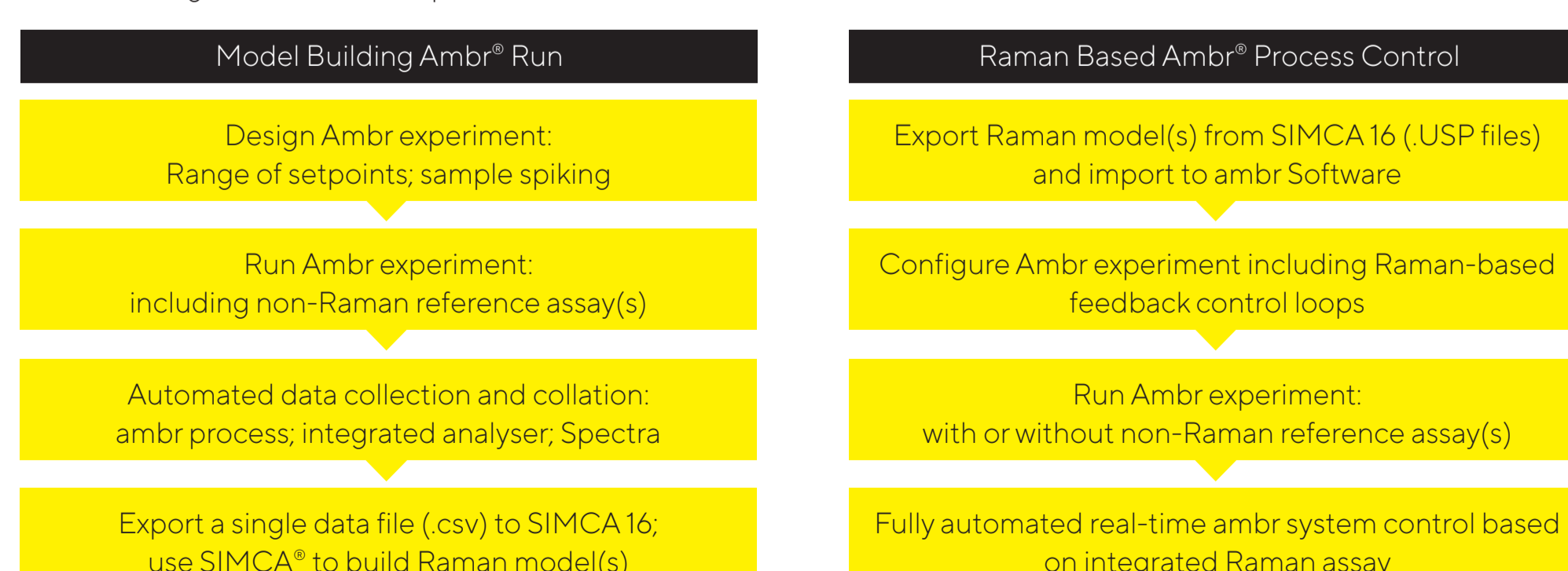

### 3. Model Building Ambr<sup>®</sup> Run

- N = 16 Ambr<sup>®</sup> 250 High Throughput bioreactors
- Sartorius' Cellca2 Process CHO producing mAb (includes glucose-containing feeds)
- Integrated Raman and Nova FLEX2 analyzers
- If daily integrated Nova FLEX2 glucose <5 g/L, automated glucose feed to 5 g/L
- >200 data points assayed by Flex2 then Raman, of which ~35 were spiked with 71 or 33.3 g/L glucose
- Data was automatically collated in the Ambr<sup>®</sup> software and one file exported by the operator
- A separate copy of SIMCA 16 software was used for model building

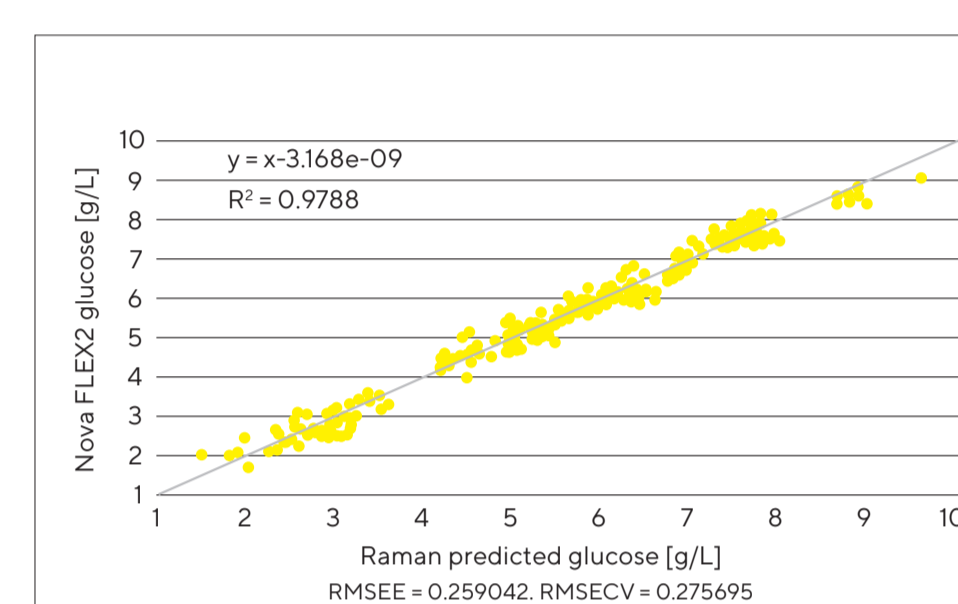


Figure 4. Glucose model excluding spiked samples (N ~ 250 total)

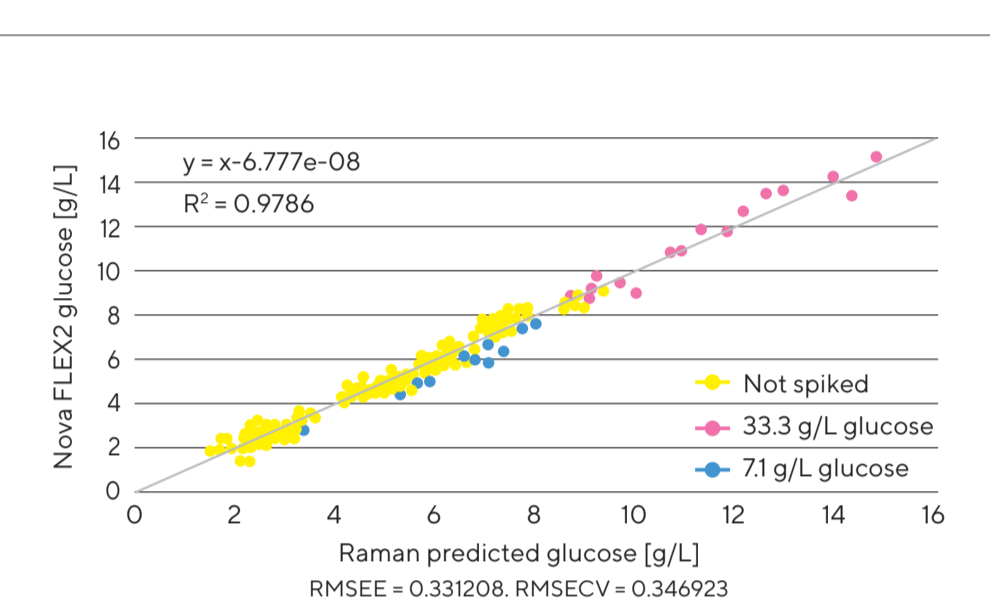


Figure 5. Glucose model, including spiked samples (N ~ 280 total)

### 4. Model Building Results

- A good Raman model was generated for glucose using only non-spiked samples (Fig. 4)
- Glucose model quality parameters are consistent with the best models in the literature
- Glucose model range was suitable for feedback control in the cell culture process investigated
- Spiking was not required to generate a good glucose model but spiking can increase model range (Fig. 5)
- Spiking has been shown to improve range and model quality for other analytes e.g. lactate (data not shown)

### 5. Raman Based Ambr<sup>®</sup> Process Control

- Prior to the run, a glucose model previously built in SIMCA<sup>®</sup> was loaded to the Ambr<sup>®</sup> software
- Sartorius' Cellca2 Process CHO producing mAb
- Integrated Raman and Nova FLEX2 analyzers
- If daily integrated Raman or FLEX2 glucose <5 g/L, automated glucose feed to 5 g/L
- N = 4 bioreactors were automatically fed glucose based on integrated Raman predictions
- N = 4 bioreactors were automatically fed glucose based on integrated Nova FLEX2 analysis

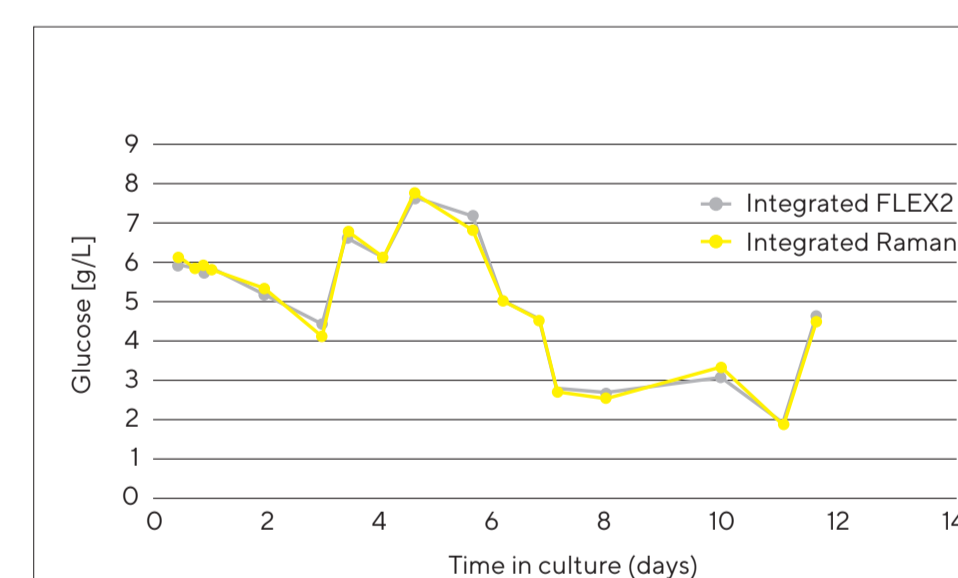


Figure 6. Glucose concentrations in a fed-batch culture, measured by both integrated FLEX2 and Raman prediction. N=1 bioreactor shown for clarity.

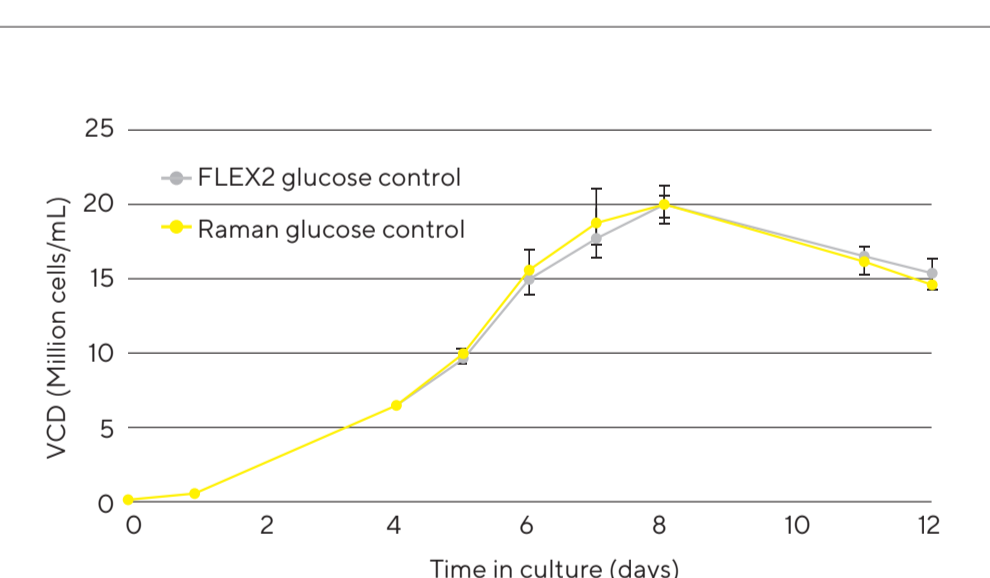


Figure 7. Cell culture VCD profiles. Glucose feeding controlled by either integrated Raman predictions (N=4) or Integrated FLEX2 assay (N=4).

### 6. Glucose Control Results

- Glucose was maintained in culture in the range of 1-8 g/L (glucose readings >5 g/L were due to glucose in feeds) (Fig. 6)
- Integrated Raman and Nova FLEX2 glucose assays were in very close agreement throughout (Fig. 6)
- Cell culture profiles (Fig. 7) were very similar for glucose control based on Raman or Nova FLEX2 assays
- BioPAT<sup>®</sup> Spectro in Ambr<sup>®</sup> enables fully automated glucose control capabilities similar to other analytical techniques (Fig. 7)
- Further work will include confirmation of the Ambr glucose model performance in Sartorius' Biostat<sup>®</sup> STR bioreactors (50-2000 L scale)

### 7. Conclusion

- Automated integration of Raman spectroscopy to Ambr<sup>®</sup> is now possible
- Fully automated data acquisition and alignment saves a significant amount of user time
- A high quality glucose model was generated covering a wide concentration range (up to 15 g/L)
- Ambr<sup>®</sup> systems enable robust model building due to a range of setpoints and spiking of samples
- Cell cultures controlled by integrated Raman or Nova FLEX2 analyzers showed very comparable glucose and VCD profiles